

INTRODUCTION


The AQ 311S Single Pan balance is of unequal arm type with 4 beams. The maximum capacity is 311 g and the readability is 0.01g. The balance comes with a stainless steel weighing pan and includes a built in platform to aid density determination by weighing samples in air and in water.

SPECIFICATIONS

Capacity	311 g
Readability	0.01g
Beam Calibration:	Front Beam 1 x 0.01g 2 nd Beam 10 x 1g 3 rd Beam 100 x 10g 4 th Beam 200 x 100 g
Damping	Magnetic
Pan Size	100mm (4") diameter Stainless Steel
Density Platform	75 mm (3") diameter Adjustable height
Overall Size (w x h x d)	370 x 300 x 137mm (14.5" x 11.8" x 5.4")
Gross Weight	2 kg. (4.4Lb)

SETTING UP

Unpack carefully. The balance consists of the main body, the pan support yoke and the stainless steel pan. Place on a firm, level surface away from fans or open windows. Position the density platform against the table. Slide all the weights to the left. Make sure the 3 largest weights fit into the notch on their beams. The front weight should point to zero. Hang the pan support yoke from the top position of the dual hook. Place the stainless steel pan on the pan support. The pointer on the right end of the beam should align with the zero mark. Adjust the thumb screw on the left end of the beam if necessary.

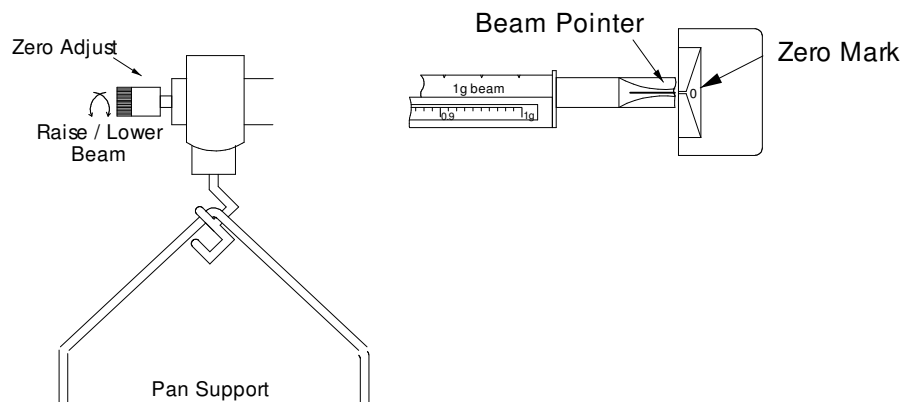

@Adam Equipment Company 2004

WEIGHING

To weigh using the AQ 311S, simply place the item to be weighed on the stainless steel pan. The beam will move up. Adjust the weights on the 4 beams until the pointer is aligned with the zero mark again. Start with the largest weight and work towards the smallest. Each weight should be moved until the beam falls below zero and moved back one notch, then the next weight is moved. The last weight moved will be the sliding weight on the front beam. Slide this weight until the beam pointer aligns with the zero mark. Read the weight by adding the values for each weight, e.g., 100 +20+5+0.27 = 125.27g.

USING A CONTAINER FOR THE SAMPLE

If an empty beaker is placed on the scale you must first determine the weight of the empty beaker. Then add the material to the beaker and weigh again. The weight of the material within the beaker is the total weight minus the weight of the beaker when empty.

FILLING TO A KNOWN WEIGHT

Place the empty beaker on the balance and determine the weight. Add the desired weight to this value to determine the final weight. Set the weights on the beam to represent this final value. Pour the material into the beaker slowly. As the desired weight is approached the beam will rise. Stop filling when the beam pointer is aligned with the zero mark.


DENSITY MEASUREMENTS

To measure the density of a solid it is necessary to find its mass in air and then in a fluid of known density. The density of the sample is then computed from:

$$\rho_s = \rho_{fl} \times \frac{m_{air}}{m_{air} - m_{fl}} \quad \rho_{fl} = \text{density of fluid}$$

If the liquid is water then ρ_{fl} is approximately 1g/cm³. See reference books to learn more about getting the best accuracy and other methods that can be used.

On the AQ 311S, this test is done by removing the pan support, raising the density platform to clear the pan and replacing the pan support. A beaker with water can be placed on the platform. Fit a piece of thread to the lower hook and adjust the weights to balance the beam. Deduct this weight from any other readings. Weigh the sample to be tested in air, either by placing it on the pan below the beaker support or by hanging from the thread on the lower hook, without the beaker in place. Weigh the sample suspended in water. Determine density from the equation above.

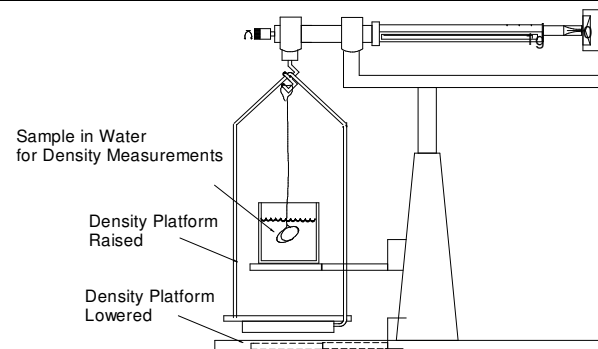

ADAM EQUIPMENT is an ISO 9001:2002 certified global organisation with more than 30 years experience in the production and sale of electronic weighing equipments. Products are sold through a world wide distributor network -supported from our company locations in the UK, USA and SOUTH AFRICA. The company and their distributors offer a full range of Technical Services such as on site and workshop repair, preventative maintenance and calibration facilities.

ADAM's products are predominantly designed for the Laboratory, Educational, Medical and Industrial Segments. The product range can be classified as follows:

- Analytical and Precision Laboratory Balances
- Top Loading Balances for Educational establishments
- Counting Scales for Industrial and Warehouse applications
- Digital Weighing/Check-weighing Scales
- High performance Platform Scales with extensive software features including parts counting, percent weighing etc.
- Digital Electronic Scales for Medical use
- Retail Scales for price computing

<p>Adam Equipment Co. Ltd. Bond Avenue Milton Keynes MK1 1SW UK</p> <p>Phone: +44 (0)1908 274545 Fax: +44 (0)1908 641339</p> <p>e-mail: sales@adamequipment.co.uk</p>	<p>Adam Equipment Inc. 26, Commerce Drive Danbury, CT 06810 USA</p> <p>Phone: +1 203 790 4774 Fax: +1 203 792 3406</p> <p>e-mail: sales@adamequipment.com</p>	<p>Adam Equipment S.A. (Pty) Ltd. P.O. Box 1422 Kempton Park 1620 Johannesburg Republic of South Africa</p> <p>Phone +27 (0)11 974 9745 Fax: +27 (0)11 392 2587</p> <p>e-mail: sales@adamequipment.co.za</p>
--	--	---

© Copyright by Adam Equipment Co. Ltd. All rights reserved. No part of this publication may be reprinted or translated in any form or by any means without the prior permission of Adam Equipment.

Adam Equipment reserves the right to make changes to the technology, features, specifications and design of the equipment without notice.

All information contained within this publication was to the best of our knowledge timely, complete and accurate when issued. However, we are not responsible for misimpressions which may result form the reading of this material.

The latest version of this publication can be found on our Website.

Visit us at www.adamequipment.com

@Adam Equipment Company 2004

AE ADAM

Adam Equipment

AQ 311S
Quadruple Beam Balances

Operator's Manual

ADAM EQUIPMENT CO. LTD.
p.n. 3866, Rev. A4, January 2004