

## **CB BALANCE**

### **In-use cover**

To attach the plastic in-use cover you must first remove the stainless steel top pan. Twist the plastic sub pan  $\frac{1}{4}$  turn to enable you to remove it from the scale. Place the in-use cover over the CB and replace the plastic sub pan securing by a  $\frac{1}{4}$  of a turn in the opposite direction, place the stainless steel top pan back on the scale. Please make sure that you do not trap the in-use cover between the CB Balance and the top pan, as this could affect the weighing accuracy.

June 2017